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The colour coding allows easy identification of the innervation of the muscles.


The peripheral nerves are demonstrated on the left hand side of the model with colour coding identifying the area of innervation of the cords of the Brachial Plexus in the upper extremity and of certain groups of the Lumbo-Sacral Plexus in the lower extremity as well as that arising from the peripheral nerves of each plexus.

The colour coding also shows the position of the respective muscle groups in the extremities and allows systematic comparison between upper and lower extremities.

The radicular or spinal nerve motor group is demonstrated on the right hand side of the model with colour coding identifying the area of innervation of certain spinal nerve roots or spinal nerves, i.e. the segment indicating muscle.


As a general rule it is a matter of one single muscle but occasionally also of various muscles and, moreover, of muscles within a number of different muscle groups of the same extremity.

The colour coding will enable a deficiency in a certain muscle within a muscle group to be attributed to a lesion in a specific motor nerve root or a specific spinal nerve.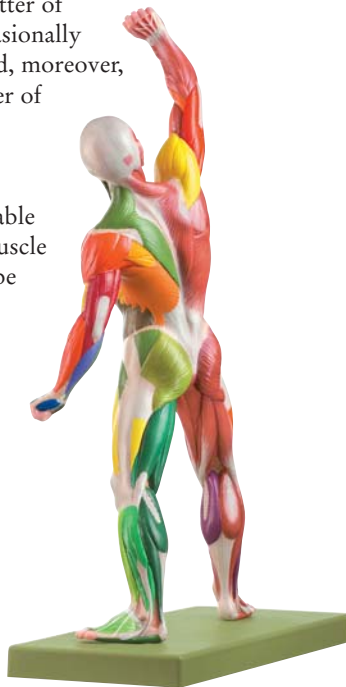


Finally the colour coding is of great assistance in allowing a differential diagnosis to be made between the involvement of the radicular or spinal nerve and the peripheral nerve in any deficiency in innervation of the musculature of an extremity.

As a result, these many teaching possibilities make the model particularly useful for various areas of applications, theoretical as well as practical from the pre-clinical stage teaching of anatomy and clinical training of orthopaedics and neurology to allied medical professions such as Physiotherapists, Sports Therapists and Masseurs.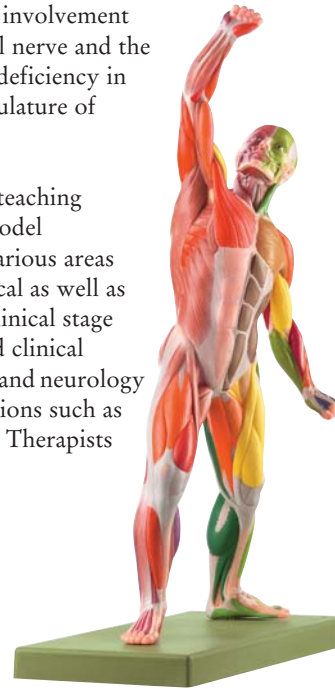

Further muscle figures from the SOMSO ANATOMY range: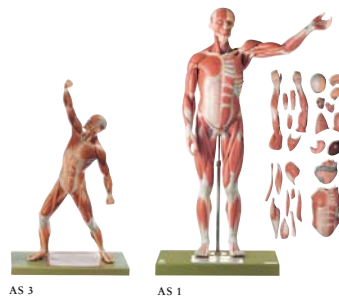

AS 1 ·
MALE MUSCLE FIGURE
about half natural
size in SOMSO-Plast
dismantleable into 27 parts.

AS 3 ·
MALE MUSCLE FIGURE
approx. quarter natural
size in SOMSO-Plast.
Not dismantleable.

Marcus Sommer
SOMSO MODELLE
Friedrich-Rückert-Str. 54
D-96450 Coburg


e-Mail: somso@somso.de
Internet: www.somso.de
Telefon (+49) (0) 95 61 / 85 74-0
Telefax (+49) (0) 95 61 / 85 74-11

Flyer A 185, © Copyright by Marcus Sommer SOMSO Modelle, August 2005

SOMSO
MODELLE
SINCE 1876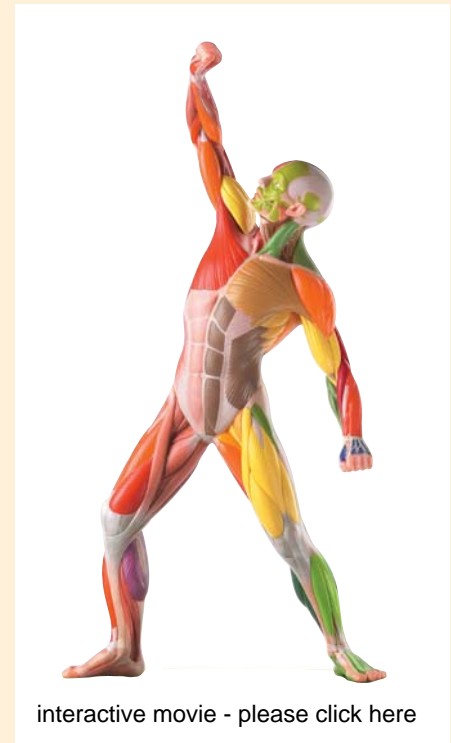

Male Muscle Figure

with colour coding for
the identification
of motor innervation

after Dr. Hans Schade

AS 3 AP/NR ·

MALE MUSCLE FIGURE WITH COLOUR CODING FOR THE IDENTIFICATION OF MOTOR INNERVATION

after Dr. Hans Schade. This model is in one piece, approx. one quarter of natural size in SOMSO-Plast. It is mounted on a plinth, removable. Height 53cm (figure 50cm) width 33cm, depth 15cm and weight 1.5 kg.

The model is as an AS3 which shows superficial musculature but with the respective muscles and muscle groups colour coded for easy identification.

Each side of the body is color coded differently.

On the left hand side of the body the muscles are colour coded corresponding to their innervation through peripheral nerves (peripheral-nerval innervation). The upper extremity colour coding is in accordance with the colour coding of FRICK-LEONHARDT-STARCK and the lower extremity has been appropriately coded using an analogous system.

On the right hand side of the body specific muscles of the upper and lower extremities the so-called "segment indicating muscles" are colour coded according to their innervation through spinal nerves roots or spinal nerves (radicular or spinal-nerval innervation). Here in accordance with the colour coding of T. von LANZ and WACHSMUTH for the spinal nerves C5-Th1 or L2-S3 respectively.

NATURE IS OUR MODEL!

SOMSO
MODELLE
SINCE 1876